



GeneAb p16 [IHC016] vs. GeneAb p16 [IHC116] vs. Ventana p16 [E6H4]

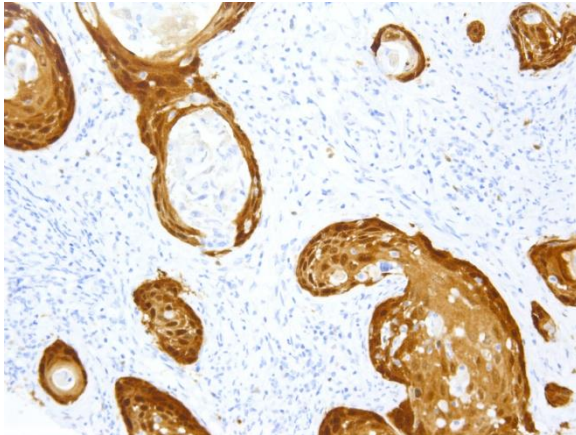


IHC016



IHC116

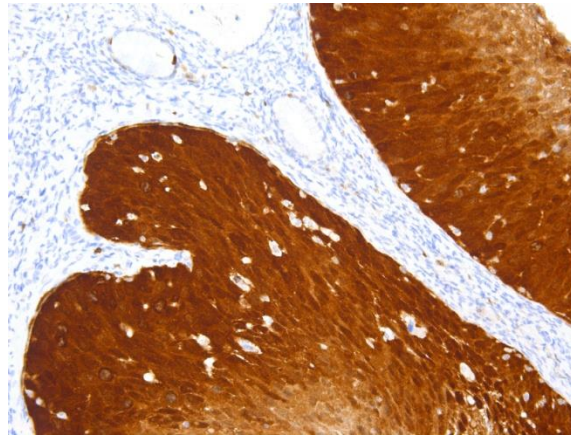

E6H4

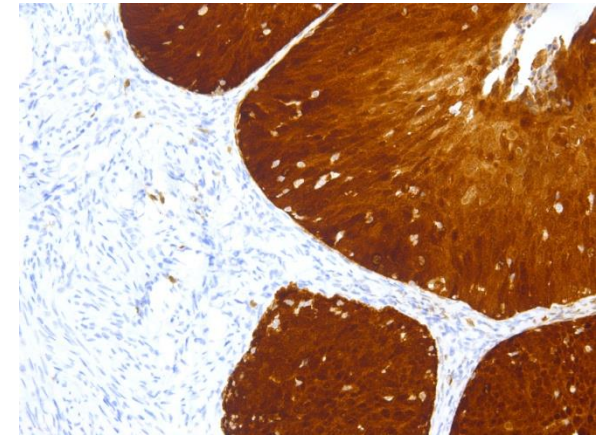

Tissue: Cervical Cancer

Staining Protocol:

Bake slides for 30 mins at 60 degrees

Dewax using dewax solution for 30s

Antigen Retrieval: HIER pretreatment using pH 9 for 30 mins

Primary Antibody Incubation: 30 mins at ambient

Post Primary Antibody Incubation: 8 mins at ambient

Polymer incubation: 8 mins at ambient

Peroxide Block: 5 mins at ambient

DAB: 10 mins at ambient

Hematoxylin: 5 mins at ambient